

## Colorimetric Microplate Assay for Total Antioxidant Power

For Research Use Only

---

### INTRODUCTION

---

Oxidative stress has been implicated in a number of diseases such as atherosclerosis<sup>1</sup>, chronic inflammatory disease<sup>2</sup>, chronic renal failure<sup>3</sup>, and cancer<sup>4</sup>. It is a condition where an imbalance exists between the production of reactive oxidizing species and the body's ability to neutralize these intermediates, resulting in cellular damage. The body has designed several physiological responses to oxidative stress including counterbalances<sup>5</sup> such as enzymes and variously functionalized molecules (see examples below) that effectively neutralize these damaging species. These antioxidants can be either water or lipid soluble, and are localized transiently throughout various tissues, cells and cell types.

Classification	Examples
Enzymes:	Superoxide Dismutase Catalase Glutathione Peroxidase
Large Molecules:	Albumin Ferritin Ceruloplasmin
Small Molecules:	Ascorbic Acid $\alpha$ -Tocopherol $\beta$ -Carotene Uric Acid

Given the multiplicity of antioxidant pathways, their centrality in the prevention of oxidant stress, and the influences of lifestyle and nutritional supplements on an individual's antioxidant capacity, it is important to be able to quantitatively measure the total antioxidant capacity or antioxidant power with biological specimens<sup>6-11</sup>.


---

### PRINCIPLES OF PROCEDURE

---

The reduction potential of the sample or standard effectively converts  $\text{Cu}^{+2}$  to  $\text{Cu}^{+1}$ , thus changing the ion's absorption characteristics. This reduced form of copper will selectively form a stable 2:1 complex with the chromogenic reagent with an absorption maximum at ca. 450 nm<sup>12-16</sup>. A known concentration of uric acid is used to create a calibration curve, with the data being expressed as mM uric acid equivalents or in  $\mu\text{M}$  copper reducing equivalents.

A study was performed to evaluate the correlation between the concentration of principal antioxidants in human serum and the value obtained for total antioxidant power using this method with the results presented in **Figure 1**. Multivariate analysis of the obtained results yields a very high significance. The results obtained with the Total Antioxidant Power Assay for a series of serum samples were also compared to the resistance to oxidation of the serum lipids in these samples. The results of this study also show that these two parameters are highly correlated. The higher the total antioxidant power, the more protected the serum lipids are from oxidation.

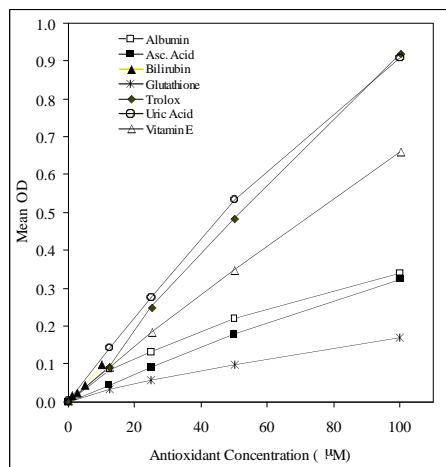

**Figure 1:** Relationship between antioxidant type and reduction from  $\text{Cu}^{+2}$  to  $\text{Cu}^{+1}$ .

### MATERIALS PROVIDED

Component	Description	Volume	Storage	Cat. No.
<b>Dilution Buffer</b>	Buffer used to dilute the standards and samples.	60 mL	4°C	TA02a
<b>Copper Solution</b>	Solution used to produce the color reaction.	5 mL	4°C	TA02b
<b>Stop Solution</b>	Solution used to stop the color reaction.	5 mL	4°C	TA02c
<b>Uric Acid Standard</b>	Lyophilized vial of 2 mM Uric Acid Standard.	1 vial	4°C	TA02d

### MATERIALS NEEDED BUT NOT PROVIDED

1. Microplate reader with a 450 nm filter
2. Adjustable micropipettes (10 – 1000 µL) and tips
3. Deionized water
4. Vortex

### STORAGE

1. Store the components of this kit at the temperatures specified on the labels.
2. Unopened reagents are stable until the indicated kit expiration date.
3. Aliquot the Uric Acid Standard after reconstitution and freeze at -80°C.

### WARNINGS AND PRECAUTIONS

1. Use aseptic technique when opening and dispensing reagents.
2. This kit is designed to work properly as provided and instructed. Additions, deletions or substitutions to the procedure or reagents are not recommended, as they may be detrimental to the assay.

### PROCEDURAL NOTES

1. Allow Dilution Buffer, Copper Solution and Stop Solution to equilibrate to room temperature for 30 minutes prior to running the assay.
2. To minimize errors in absorbance measurements due to handling, wipe the exterior bottom of the microplate wells with a lint-free paper towel prior to inserting into the plate reader.

---

## SAMPLE PREPARATION

---

All samples should be stored at  $-70^{\circ}\text{C}$  *at all times* prior to assay. Apply the samples to the assay immediately upon thawing. Samples with a uric acid equivalent concentration greater than 2.0 mM should be diluted with pH 7.0 PBS prior to assay. Sample preparation is subject to the discretion and approval of the principal investigator, and changes may occur to optimize or better complement experimental design.

**NOTE:** Ethylene diamine tetraacetic acid (EDTA) is detrimental to the function of the assay and *must not be present at any concentration* in any sample type.

**Urine:** Collected samples may be assayed directly or diluted with pH 7.0 PBS where appropriate.

**Tissue Lysate:** Homogenize tissue sample on ice with ice-cold pH 7.0 PBS then centrifuge at  $3000 \times g$  for 12 minutes at  $4^{\circ}\text{C}$ . Aliquot the supernatant for storage at  $-70^{\circ}\text{C}$ , protein determination and subsequent assays.

**Cell Culture:** Wash cells 2-3 times with ice-cold pH 7.0 PBS prior to lysis. Lyse cells by homogenization or sonication with ice-cold pH 7.0 PBS then centrifuge at  $3000 \times g$  for 12 minutes at  $4^{\circ}\text{C}$ . Aliquot the supernatant for storage at  $-70^{\circ}\text{C}$ , protein determination and subsequent assays.

**Plasma:** Collect blood with sodium citrate and centrifuge at  $3000 \times g$  for 12 minutes at  $4^{\circ}\text{C}$ . Remove the plasma and aliquot for storage at  $-70^{\circ}\text{C}$ , protein determination and subsequent assays. Sample results can alternatively be expressed in mL of plasma.

**Food:** This can vary considerably from one food to the next and requires foresight and discretion from the principal investigator. In many cases liquid food samples such as juice and tea can be assayed directly without any processing. Liquid samples that contain high amounts of protein and/or fiber or solid foods should be processed as indicated below.

Solid Food: Homogenize food sample on ice in ice-cold pH 7.0 PBS then centrifuge at  $3000 \times g$  for 12 minutes at  $4^{\circ}\text{C}$ . Aliquot the supernatant for storage at  $-70^{\circ}\text{C}$ , protein determination and subsequent assays.

Liquid Food: Centrifuge at  $3000 \times g$  for 12 minutes at  $4^{\circ}\text{C}$ . Aliquot the supernatant for storage at  $-70^{\circ}\text{C}$ , protein determination and subsequent assays.

---

## STANDARD CURVE PREPARATION

---

**PLEASE NOTE:** The Uric Acid Standard vial is under vacuum.

### TO RECONSTITUTE THE URIC ACID STANDARD:

**Preferred Method** – Add 1.5 mL of deionized water directly to the Uric Acid Standard vial with a needle and syringe by puncturing the rubber stopper with the needle. Vortex the vial for 30-60 seconds, or until the standard is completely dissolved. The concentration is now 2.0 mM. Slowly remove the rubber stopper, allowing the vacuum seal to break and proceed to **Table 1** for the standard curve dilutions.

**Alternative Method** – Not Recommended - WARNING!!! The solid uric acid can become airborne due to the opening process or due to the turbulence created during pipetting and reconstitution. Loss of material will greatly affect the accuracy of your assay – this method is performed at the user's risk and discretion. Slowly lift a corner of the rubber stopper until the vacuum seal is broken. Now remove the stopper and slowly add 1.5 mL of DI  $\text{H}_2\text{O}$  to the vial. Recap the vial with the same rubber stopper and Parafilm or tape the stopper to the vial then vortex for 30-60 seconds, or until the standard is completely dissolved. The concentration of this standard is now 2.0 mM. Carefully remove the stopper and proceed to **Table 1** for the standard curve dilutions.

The reconstituted standard can be aliquoted and stored at  $-70^{\circ}\text{C}$  for up to one year.

Use the following table to dilute the 2 mM Uric Acid Standard and construct a six-point standard curve.

**Table 1:** Standard Curve Preparation

Standard	Uric Acid Conc. (mM)	Vol. of Deionized Water ( $\mu$ L)	Transfer Volume ( $\mu$ L)	Transfer Source	Final Volume ( $\mu$ L)
S <sub>5</sub>	2.0	1500	-	2 mM Stock	1000
S <sub>4</sub>	1.0	500	500	S <sub>5</sub>	500
S <sub>3</sub>	0.5	500	500	S <sub>4</sub>	500
S <sub>2</sub>	0.25	500	500	S <sub>3</sub>	500
S <sub>1</sub>	0.125	500	500	S <sub>2</sub>	1000
S <sub>0</sub>	0	500	-	-	500

---

**ASSAY PROCEDURE**


---

1. Dilute both Samples and Standards 1:40 in the provided Dilution Buffer (e.g. 15  $\mu$ L serum + 585  $\mu$ L Dilution Buffer). See **Scheme I** for a suggested plate layout.
2. Place 200  $\mu$ L of diluted Samples or Standards in each well. Reagent Blanks (BLK) should contain Dilution Buffer in place of Standard or Sample.
3. Read the plate at 450 nm for a reference measurement.
4. Add 50  $\mu$ L of Copper Solution to each well and incubate for 3 minutes at room temperature.
5. Add 50  $\mu$ L of Stop Solution.
6. Read the plate a second time at 450 nm.

**Scheme I:** Sample Plate Layout

	1	2	3	4	5	6	7	8	9	10	11	12
<b>A</b>	S <sub>0</sub>	S <sub>0</sub>	U <sub>3</sub>	U <sub>3</sub>	U <sub>11</sub>	U <sub>11</sub>	U <sub>19</sub>	U <sub>19</sub>	U <sub>27</sub>	U <sub>27</sub>	U <sub>35</sub>	U <sub>35</sub>
<b>B</b>	S <sub>1</sub>	S <sub>1</sub>	U <sub>4</sub>	U <sub>4</sub>	U <sub>12</sub>	U <sub>12</sub>	U <sub>20</sub>	U <sub>20</sub>	U <sub>28</sub>	U <sub>28</sub>	U <sub>36</sub>	U <sub>36</sub>
<b>C</b>	S <sub>2</sub>	S <sub>2</sub>	U <sub>5</sub>	U <sub>5</sub>	U <sub>13</sub>	U <sub>13</sub>	U <sub>21</sub>	U <sub>21</sub>	U <sub>29</sub>	U <sub>29</sub>	U <sub>37</sub>	U <sub>37</sub>
<b>D</b>	S <sub>3</sub>	S <sub>3</sub>	U <sub>6</sub>	U <sub>6</sub>	U <sub>14</sub>	U <sub>14</sub>	U <sub>22</sub>	U <sub>22</sub>	U <sub>30</sub>	U <sub>30</sub>	U <sub>38</sub>	U <sub>38</sub>
<b>E</b>	S <sub>4</sub>	S <sub>4</sub>	U <sub>7</sub>	U <sub>7</sub>	U <sub>15</sub>	U <sub>15</sub>	U <sub>23</sub>	U <sub>23</sub>	U <sub>31</sub>	U <sub>31</sub>	U <sub>39</sub>	U <sub>39</sub>
<b>F</b>	S <sub>5</sub>	S <sub>5</sub>	U <sub>8</sub>	U <sub>8</sub>	U <sub>16</sub>	U <sub>16</sub>	U <sub>24</sub>	U <sub>24</sub>	U <sub>32</sub>	U <sub>32</sub>	U <sub>40</sub>	U <sub>40</sub>
<b>G</b>	U <sub>1</sub>	U <sub>1</sub>	U <sub>9</sub>	U <sub>9</sub>	U <sub>17</sub>	U <sub>17</sub>	U <sub>25</sub>	U <sub>25</sub>	U <sub>33</sub>	U <sub>33</sub>	U <sub>41</sub>	U <sub>41</sub>
<b>H</b>	U <sub>2</sub>	U <sub>2</sub>	U <sub>10</sub>	U <sub>10</sub>	U <sub>18</sub>	U <sub>18</sub>	U <sub>26</sub>	U <sub>26</sub>	U <sub>34</sub>	U <sub>34</sub>	BLK	BLK

---

**CALCULATIONS**


---

1. Calculate the net absorbance by subtracting the absorbance readings from step 3 from those in step 6.
2. Plot the net absorbance vs. uric acid concentration. A typical standard curve is shown in **Figure 2**.

Samples can be expressed in terms of  $\mu$ M uric acid equivalents by solving for “x” from your generated standard curve equation (see example below). To express values in  $\mu$ M copper reducing equivalents (CRE’s), multiply the “x” value coefficient by 2189  $\mu$ M.

$$y = mx$$


where: y = y-axis value (OD readings)

x = x-axis value (Uric Acid concentration)

m = slope

Given from the standard curve provided on the next page:  $m = 0.2264$

If “y” has an OD value of 0.23 then “x” equals 1.0 mM uric acid equivalents. To express as CRE’s, multiply the mM equivalents of uric acid by 2,189 to obtain the CRE value.

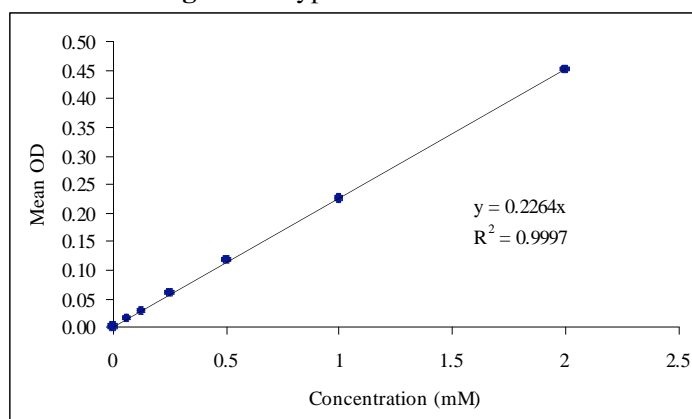**Figure 1: Typical Standard Curve**

---

**REFERENCES**

---

1. Brook, R.; (2008) *Clin. Sci. (Lond)*. **115**:175-187
2. Cerutti, P. and Trump, B.; (1991) *Cancer Cell* **3**:1-7
3. Cottone, S., et al.; (2008) *J. Nephrol.* **21**:175-179
4. Trachootham, D., et al.; (2008) *Antioxid. Redox Signal.* **10**:1343-1374
5. Frei, B., et al.; (1992) *Molecular Biology of Free Radical Scavenging System* 23-45
6. Ciencewicki, J., et al.; (2008) *J. Allergy Clin. Immunol.* **122**:456-468
7. Kieth, M., et al.; (1998) *J. Am. Coll. Cardio.* **31**:1352-1362
8. Van-Zoeren-Grobben, et al.; (1997) *Acta Paediatrica* **86**:1356-1362
9. Allard, J.P., et al.; (1998) *Am. J. Clin. Nutr.* **67**:143-147
10. Santini S.A., et al.; (1997) *Diabetes* **46**:1853-1858
11. Fatokun, A., Stone, T. and Smith, R.; (2008) *Front. Biosci.* **13**:3288-3311
12. Apak, R., et al.; (2004) *J Agric Food Chem* **52**:7970-7981
13. Apak, R., et al.; (2005) *Free Radic Res* **39**:949-961
14. Apak, R., et al.; (2007) *Molecules* **12**:1496-1547
15. Apak, R., et al.; (2008) *Microchim Acta* **160**:413-419
16. Apak, R., et al.; (2008) *Anal. Chim. Acta* **630**:28-39

---

**DISCLAIMER**

---

This information is believed to be correct but does not purport to be all-inclusive and shall be used only as a guide. Oxford Biomedical Research, Inc. shall not be held liable for any damage resulting from handling or from contact with the above product. See catalog for additional terms and conditions of sale.

---

**ORDERING INFORMATION**

---

For additional kits or a complete catalog please call 800-692-4633.

---

**TECHNICAL SUPPORT**

---

If you need technical information or assistance with assay procedures, call our Technical Support Department at 800-692-4633 or 248-852-8815. Our staff will be happy to answer your questions about this or any other product in the Oxford Biomedical line.

---

**GUARANTEE AND LIMITATION OF REMEDY**

---

Oxford Biomedical Research, Inc. makes no guarantee of any kind, expressed or implied, which extends beyond the description of the materials in this kit, except that these materials and this kit will meet our specifications at the time of delivery. Buyer's remedy and Oxford Biomedical Research, Inc.'s sole liability hereunder is limited to, at Oxford Biomedical Research, Inc.'s option, refund of the purchase price of, or the replacement of, material that does not

meet our specification. By acceptance of our products, Buyer indemnifies and holds Oxford Biomedical Research, Inc. harmless against, assumes all liability for the consequence of its use or misuse by the Buyer, its employees, or others. Said refund or replacement is conditioned on Buyer notifying Oxford Biomedical Research, Inc. within thirty (30) days of the receipt of product. Failure of Buyer to give said notice within thirty (30) days of receipt of product shall constitute a waiver by the Buyer of all claims hereunder with respect to said material(s).

Oxford Biomedical Research, Inc.  
P.O. Box 522  
Oxford, MI 48371 U.S.A.

Orders: 800-692-4633  
Technical Service: 248-852-8815  
Fax: 248-852-4466  
E-mail: [info@oxfordbiomed.com](mailto:info@oxfordbiomed.com)

Made in the U.S.A.