

Fluorometric Microplate Assay for 2-Thiobarbituric Acid Reactive Substances (TBARS)

For Research Use Only

INTRODUCTION

2-ThioBarbituric Acid Reactive Substances (TBARS) are naturally present in biological specimens and include lipid hydroperoxides and aldehydes which increase in concentration as a response to oxidative stress.^{1,2} TBARS assay values are usually reported in malonaldehyde (malondialdehyde, MDA) equivalents, a compound that results from the decomposition of polyunsaturated fatty acid lipid peroxides. The TBARS assay is a well-recognized, established method for quantifying these lipid peroxides, although it has been criticized for its reactivity towards other compounds other than MDA.³ This kit contains modifications from the traditional TBARS assay to reduce background and the generation of TBA reactive substances to give a straightforward, reproducible and consistent method for analyzing biological samples for lipid peroxidation products.

PRINCIPLES OF PROCEDURE

This assay is based on the reaction of a chromogenic reagent, 2-thiobarbituric acid, with MDA at 25°C. One molecule of MDA reacts with 2 molecules of 2-thiobarbituric acid *via* a Knoevenagel-type condensation to yield a chromophore with absorbance maximum at 532 nm, as shown below in Figure 1:

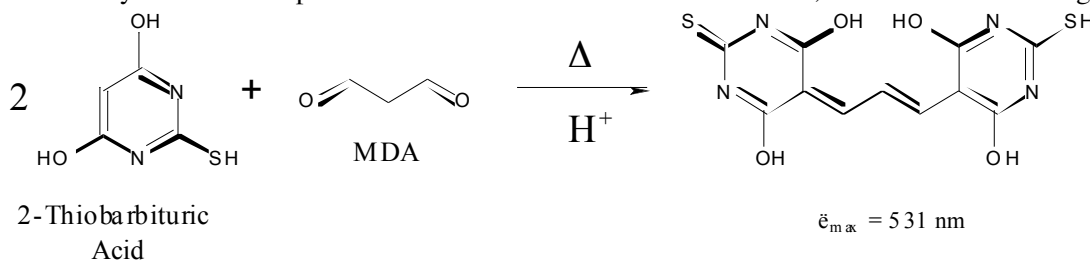


Figure 1: Reaction between 2-thiobarbituric acid and MDA under acidic conditions.

MATERIALS PROVIDED

Component	Description	Amount	Storage	Cat. No.
Indicator	2-Thiobarbituric Acid	2 x 0.5 g	4°C	FR45a
Acid Reagent	10% Acid Solution in Dimethylsulfoxide	40 mL	4°C	FR45b
MDA Standard	10 mM Malonaldehyde Tetrabutylammonium Salt	100 μ L	4°C	FR45c
Microplate	96-well Microplate	2 plates	4°C	FR45d

MATERIALS NEEDED BUT NOT PROVIDED

1. Fluorometric plate reader capable of exciting the sample at 532 nm and reading at 585 nm.
2. Polypropylene microcentrifuge tubes
3. Deionized Water (dH_2O)
4. Adjustable micropipettes (10 – 1,000 μ L) and tips

STORAGE

1. The reagents are stable until the indicated kit expiration date if handled and stored properly.
2. When not in use, store the kit at 4°C for up to one year.
3. MDA standards should be used within 24 hours of preparation.
4. The Indicator Solution (combined 2-TBA and Acid Solution) can be stored at 4°C for one week.

WARNINGS AND PRECAUTIONS

1. Use aseptic techniques when opening and dispensing reagents.
2. Wear gloves and safety glasses when performing this assay, as the acid used is corrosive.
3. In case of accidental exposure to 2-TBA or Acid Reagent, thoroughly wash the exposed area with soap and water.
4. This kit is designed to work properly as provided and instructed. Additions, deletions or substitutions to the procedure or reagents are not recommended, as they may be detrimental to the assay.

PROCEDURAL NOTES

1. Do not leave the reagent bottles open. Replace the caps as soon as the desired volume is removed.
2. To minimize error due to handling, wipe the exterior bottom of the microplate wells with a lint-free paper towel.
3. **Sample Blanks:** It is HIGHLY recommended that, for each sample analyzed, a sample blank is also performed to account for background interferences. While Oxford's TBARS method has relatively low background, some background interferences are inherent in biological samples and cannot be avoided. The sample blank should contain the sample plus the Acid Reagent alone without TBA indicator. Following this procedure will result in more accurate, precise and reliable TBARS measurements.
4. There are sufficient standards and reagents to test 40 samples and 40 sample blanks in duplicate. Samples and sample blanks should be run together on the same plate using the same standard curve.

REAGENT PREPARATION

1. **Acid Reagent:** This solution may need to be thawed upon removal from 4°C. Allow to sit at room temperature for one hour. It is ready to use once thawed.
2. **Indicator Solution:** Add 10 mL of the Acid Reagent to the powdered contents of one vial of Indicator and shake until completely dissolved. One vial is sufficient for standards and 20 samples.
3. **20 nM MDA Standard Stock:** First, dilute the 10 mM MDA Standard 1:500 in dH₂O by adding 20 µL of 10 mM MDA to 9.98 mL dH₂O. Then dilute this solution 1:1,000 in dH₂O by adding 10 µL of 20 µM MDA to 9.99 mL dH₂O. Prepare immediately prior to use.

SAMPLE STABILITY

Studies at OBR show that this kit provides optimum results with urine that is less than 48 hours old, and is preferably run immediately after sample collection. If this is not possible, samples should be frozen at -70°C to prevent loss of MDA and HAE⁴ and sample oxidation. Samples should **not** be stored at -20°C. Samples should not be refrozen and should be protected from light to avoid photooxidation.

SAMPLE PREPARATION

When working with plasma, the sample should be deproteinated with an acid. Centrifuge and use the supernatant to perform the assay. This solution may appear cloudy after the reaction, and can be clarified by passing through a 0.2 µ syringe filter.

When working with urine, colored compounds contribute to the signal measured at 532 nm. This interference can be removed by running a sample blank with each sample.

Urine

1. Urine samples can be used directly and should be assayed immediately. If the assay is to be performed on a different day, the sample should be stored at -70°C .

Plasma and Serum

1. Collect blood samples and process immediately per the collection tube instructions.
2. Prepare a saturated solution of ammonium sulfate.
3. Add 100 μL of saturated ammonium sulfate to 0.5 mL of serum or plasma in a test tube or microcentrifuge tube.
4. Add 35 mg TCA (trichloroacetic acid) to each sample and vortex. A cloudy precipitate should form.
5. Centrifuge the tubes and transfer the supernatant to a clean tube. Plasma and serum samples can be run without dilution. Samples are now ready for assay.

STANDARD CURVE PREPARATION

Malondialdehyde is provided as a solution of the malondialdehyde tetrabutylammonium (MDA-TBA) salt in a slightly basic buffer because MDA itself is not stable. When mixed with the acidic Indicator Solution, the MDA-TBA molecule is acidified and generates MDA quantitatively.

Please see the **Reagent Preparation** section for preparing the 20 nM MDA Standard Stock.

Table 1: Fluorometric Standard Curve Preparation

Standard	MDA Conc. (nM)	Vol. of dH ₂ O (μL)	Vol. of 20 nM MDA Stock (μL)
S ₀	0	400	-
S ₁	0.5	390	10
S ₂	1.0	380	20
S ₃	2.5	350	50
S ₄	5.0	300	100
S ₅	10.0	200	200
S ₆	15.0	100	300
S ₇	20.0	-	400

ASSAY PROCEDURE

Free MDA

1. **Preparation of Standards and Samples:** Add each of the following reagents into microcentrifuge tubes and mix well.
 - **Standards:** 200 μL of standard and 200 μL of Indicator Solution.
 - **Samples:** 200 μL of sample and 200 μL of Indicator Solution.
 - **Blanks:** 200 μL of sample and 200 μL of Acid Reagent.
2. After the standards, samples and blanks have been mixed; allow them to react for 45 minutes at room temperature.
3. Transfer 150 μL of each solution to the black microplate and excite the sample at 532 nm and read at 585 nm. See Scheme 1 on the following page for a sample plate layout.

Total MDA


1. Prepare samples and standards exactly as above, but heat sample at 65°C for 45 minutes, then follow step 3 as above.

Scheme 1: Sample Plate Layout

	1	2	3	4	5	6	7	8	9	10	11	12
A	S ₀	S ₀	SPL ₁	SPL ₁	SPL ₅	SPL ₅	SPL ₉	SPL ₉	SPL ₁₃	SPL ₁₃	SPL ₁₇	SPL ₁₇
B	S ₁	S ₁	SB ₁	SB ₁	SB ₅	SB ₅	SB ₉	SB ₉	SB ₁₃	SB ₁₃	SB ₁₇	SB ₁₇
C	S ₂	S ₂	SPL ₂	SPL ₂	SPL ₆	SPL ₆	SPL ₁₀	SPL ₁₀	SPL ₁₄	SPL ₁₄	SPL ₁₈	SPL ₁₈
D	S ₃	S ₃	SB ₂	SB ₂	SB ₆	SB ₆	SB ₁₀	SB ₁₀	SB ₁₄	SB ₁₄	SB ₁₈	SB ₁₈
E	S ₄	S ₄	SPL ₃	SPL ₃	SPL ₇	SPL ₇	SPL ₁₁	SPL ₁₁	SPL ₁₅	SPL ₁₅	SPL ₁₉	SPL ₁₉
F	S ₅	S ₅	SB ₃	SB ₃	SB ₇	SB ₇	SB ₁₁	SB ₁₁	SB ₁₅	SB ₁₅	SB ₁₉	SB ₁₉
G	S ₆	S ₆	SPL ₄	SPL ₄	SPL ₈	SPL ₈	SPL ₁₂	SPL ₁₂	SPL ₁₆	SPL ₁₆	SPL ₂₀	SPL ₂₀
H	S ₇	S ₇	SB ₄	SB ₄	SB ₈	SB ₈	SB ₁₂	SB ₁₂	SB ₁₆	SB ₁₆	SB ₂₀	SB ₂₀

CALCULATIONS

1. Average the Relative Light Unit (RLU) values obtained for all duplicated wells.
2. Plot a standard curve using the RLU value for each Standard versus the MDA concentration for each Standard. The equation of the line can be found using a linear fit method.
3. Subtract the RLU for each Sample Blank from the Sample RLU to obtain a Net RLU.
4. Calculate the MDA concentration for each Sample using the Net RLU value and the equation generated by the Standard Curve. If the Samples were diluted, the result must be multiplied by the dilution factor. If you are testing serum or plasma samples and have followed the deproteinization step above, your dilution factor is 1.2.

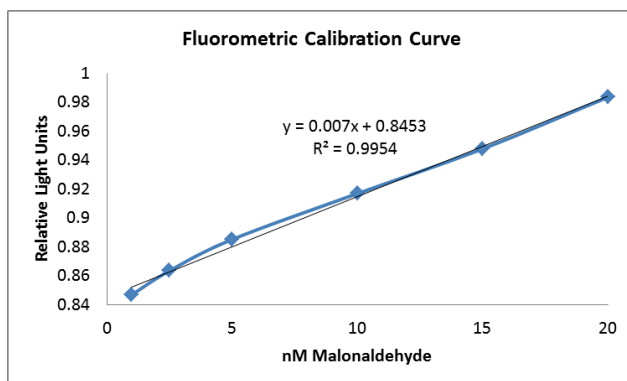Figure 2: Typical Standard Curve**LIMIT OF DETECTION**

The limit of detection for the fluorometric assay has been determined to be 5.0 nM.

PERFORMANCE LIMITATIONS

1. Although the standards in this assay will usually appear water clear, the samples may become colored. This is due to the formation of additional chromophores that absorb at various wavelengths other than 532 nm and will usually not interfere with the A_{532} signal.
2. In setting up this assay for the first time on a particular biological sample, the kinetics of color development in the samples should be followed in comparison with those of the MDA standards. The A_{532} of the standards should reach a plateau after approximately 15 minutes and then remain. If the A_{532} signal continues to increase after the standards have achieved a stable color, the researcher should be concerned that interfering non-TBARS related reactions are occurring in the sample.

REFERENCES

1. Armstrong, D.; Browne, R. (1994) *Free Rad. Diag. Med.* **366**; 43-58.
2. Botsoglou, N.A., *et al.*; (1994) *J. Agric. Food Chem.* **42**; 1931-1937.
3. Yagi, K. (1998) *Free Rad. Antiox. Prot.*, **108**; 101-106.
4. Moore, K., Roberts, L.J., (1998) *Free Rad. Res.* **28**; 659-671.
5. Liu, J., *et al.*; (1997) *Analyt. Biochem.* **245**;161-166.
6. Carbonneau, M.A., *et al.*; (1991) *Clin. Chem.* **37**; 1423-1429.
7. Mattson, J.P., *et al.*; (2002) *Pathophysiology* **8**; 215-221.
8. Calvo, J.R., *et al.*; (2001) *J. Cell Biochem.* **80**; 461-70.
9. Janciauskiene, S. and Bo Ahren; (2000) *Biochem. Biophys. Res. Com.* **267**; 619-625.
10. Montilla-Lopez, P., *et al.*; (2002) *European J. Pharmacology* **451**; 237-243.

DISCLAIMER

This information is believed to be correct but does not purport to be all-inclusive and shall be used only as a guide. Oxford Biomedical Research, Inc. shall not be held liable for any damage resulting from handling or from contact with the above product. See catalog for additional terms and conditions of sale.

ORDERING INFORMATION

For additional kits or a complete catalog please call 800-692-4633.

TECHNICAL SUPPORT

If you need technical information or assistance with assay procedures, call our Technical Support Department at 800-692-4633 or 248-852-8815. Our staff will be happy to answer your questions about this or any other product in the Oxford Biomedical line.

GUARANTEE AND LIMITATION OF REMEDY

Oxford Biomedical Research, Inc. makes no guarantee of any kind, expressed or implied, which extends beyond the description of the materials in this kit, except that these materials and this kit will meet our specifications at the time of delivery. Buyer's remedy and Oxford Biomedical Research, Inc.'s sole liability hereunder is limited to, at Oxford Biomedical Research, Inc.'s option, refund of the purchase price of, or the replacement of, material that does not meet our specification. By acceptance of our products, Buyer indemnifies and holds Oxford Biomedical Research, Inc. harmless against, assumes all liability for the consequence of its use or misuse by the Buyer, its employees, or others. Said refund or replacement is conditioned on Buyer notifying Oxford Biomedical Research, Inc. within thirty (30) days of the receipt of product. Failure of Buyer to give said notice within thirty (30) days of receipt of product shall constitute a waiver by the Buyer of all claims hereunder with respect to said material(s).

Oxford Biomedical Research, Inc.
Orders: 800-692-4633
Technical Service: 248-852-8815
E-mail: info@oxfordbiomed.com
Made in the U.S.A.